

# Bulb Replacement for Lasico model 7000 Crystalscopes with non slotted Rack Lifter (17)

1. Remove the microscope tube (8)  
After loosening clamp screw (51)  
pull the tube out



2. Remove the clamp release (5) after  
opening the pivot screw (33)



3. Remove two hex bolts (50) which fasten  
the bridge assembly (2) to the  
Crystalscope housing and remove  
the bridge assembly (2)



4. Unsolder the wire from the power jack (35)



5. Loosen the set screw of the lock ring (34) and pull the lock ring out




6. Open and remove the plug screw (21) of the lamp housing



7. Using pliers (careful not to damage the insulation) and pull the cable out while pushing it from above down through the rack lifter (17)



8. Subsequently pull the lamp housing (25) out of the Crystalscope

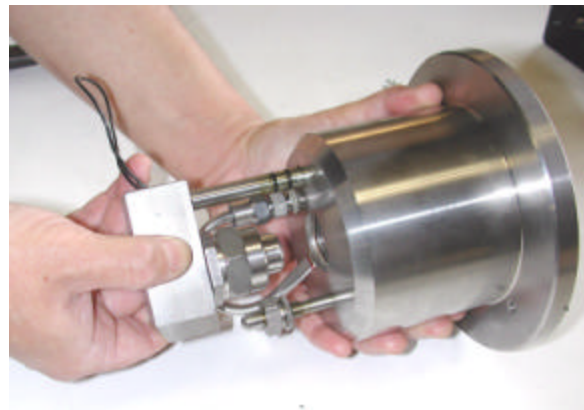

9. Unscrew the lens housing (27)



10. Replace bulb



11. Reassemble in reverse order



2451 Riverside Drive, Los Angeles, Ca.90039 -USA  
Phone: (323) 662-2128 Fax: (323) 662-0904 E-Mail : [lasico@worldnet.att.net](mailto:lasico@worldnet.att.net)  
Website: [www.lasico.com](http://www.lasico.com) [www.planimeters.com](http://www.planimeters.com)